

Sampling Instructions for "Basic" City Water Testing Kit

Carefully Read Instructions Before Collecting Your Samples

- This kit will test for Nitrate, Nitrite, Fluoride, Chloride, Sulfate, Lead, Copper, Sodium, Calcium, Magnesium, Hardness, pH, Total Dissolved Solids
- Cooler contains: Blue Gel Ice Pack(s), 2 Large Round Bottles, 1 Small Round Bottle, these Instructions, and a Laboratory Form
- Identify whether or not the required shipping arrangements can be met before collecting the samples
- One Sampling Kit to be used for One Sampling Point

1. Take out the Blue Gel Ice Pack from the cooler and freeze it for at least 8 hours prior to sample collection. Return to the cooler just prior to sampling and/or shipping.
2. Determine the sampling point (kitchen tap, bathroom tap, etc.).
3. Remove any aerators, purifiers, or other devices from the tap/spigot you are collecting samples from.
4. Run the cold water for approximately 15 minutes, then adjust the flow to a thin stream (about the diameter of a pencil).

Do not forget to write the Name of Sampling Point, Date and Time of sampling on each bottle label

5. First, obtain one large round bottle (Test for Lead, Copper) Fill the bottle to the neck and cap securely. Place the bottle in the cooler.
6. Next, obtain the second large round bottle (Test for Nitrate, Nitrite, Fluoride, Chloride, Sulfate, pH, Total Dissolved Solids). Fill the bottle to the neck and cap securely. Place the bottle in the cooler.
7. Then obtain the small round bottle (Test for Sodium, Calcium, Magnesium, Hardness). Fill the bottle to the neck and cap securely. Place the bottle in the cooler.
8. Complete the enclosed Laboratory Form, indicating the Name of Sampling Point (kitchen tap, bathroom tap, etc.), Date and Time of sample collection for each bottle. Retain a copy of the Laboratory Form for your records and return the completed form to the cooler.
10. Return the blue gel ice pack(s) to the cooler and prepare the cooler for delivery/shipment.

Shipping Arrangements

*****The collected samples MUST be received by the laboratory within 24 hours of sample collection*****

11. If you need to ship the collected samples back to the laboratory, be sure that the carrier (UPS, FedEx, etc.) will deliver the samples within 24 hours from the time of the sample collection. It is suggested that the sample collection take place as close to the time of shipping as possible to ensure the time constraint will be met.
12. If you will deliver the collected samples to the laboratory, please call us at 1-800-259-9532 to make necessary arrangements.

Laboratory address: Advanced Analytical Technologies, Inc.
 37 Ramland Road
 Orangeburg, NY 10962

Please call us at 1-800-259-9532 if you have any questions regarding sample collection or our services.